



C Y C L E X - D S A M P L I N G G U I D E

SAMPLING PROCEDURE

1. Remove blue pin from the bottom of the cycllex-d.
2. Simply connect one end of tubing (5 feet or less) to the bottom of the cycllex-d (No adapters are necessary).
3. Connect the other end of tubing to a pump precalibrated to a flow rate of 20 lpm.
4. Remove red cap from the top of the cycllex-d.
5. Turn pump on and pull a sample for 10 minutes, depending on environment (See "Sampling Intervals").
6. When sampling is complete, replace blue pin and red cap to the bottom and top of the cycllex-d.

Sampling Intervals

Clean (Outdoor/Indoor)
 High Activity (Indoor)
 Inner Wall Samples

Sampling Time

10 minutes
 5-8 minutes
 2-5 minutes

